BioClean™ GOGGLES

INSTRUCTIONS FOR USE

P-IFU-GOG-V5

Available in 29 languages
v. 2024-05

Please click on the box below to select your preferred language.

Europe, Middle East and Africa Region

Ansell Healthcare Europe NV
Riverside Business Park
Blvd International, 55
1070 Brussels, Belgium

UK Importer / Authorized Rep:

Nitritex Ltd
Ground Floor, 15 Kings Court,
Willie Snaith Road, Newmarket,
Suffolk, CB8 7SG, United Kingdom

www.ansell.com

Ansell, ® and ™ are trademarks owned by Ansell Limited or one of its affiliates. © 2024 Ansell Limited. All Rights Reserved.

CONTENTS

Please click on the content below to navigate to your selected page.

SYMBOLS & PICTOGRAMS

EN166:2001

(01)

(02)

(03)

(04)

(05)

EU Declaration of conformity available to download at - www.ansell.com/regulatory | EU Notified Body: INTERTEK Italia S.p.A. Via Guido Miglioli 2/A, 20063 Cernusco sul Naviglio - Milano (MI), Italy. Notified Body. No. 2575 | **UKCA Approved Body:** ITS Testing Services (UK) Ltd, Centre Court, Meridian Business Park, Leicester, UK. Approved Body No. AB0362.

BioClean™ goggles are intended to protect the eyes from the following hazards defined by the fields of use in EN 166: 2001

| EN 166 Field of use | BCGS1 | BVGS | BCAG | BCAP | BCAH |
|---|-------|------|------|------|------|
| 3 Liquids | ✓ | ✓ | X | ✓ | ✓ |
| 4 Large dust particles | ✓ | ✓ | X | X | X |
| B Medium impact energy | ✓ | ✓ | ✓ | ✓ | ✓ |
| BT Medium impact energy at extremes of temperatures | X | X | ✓ | ✓ | ✓ |
| N Fogging | X | X | ✓ | ✓ | ✓ |
| 1 Optical class 1 | ✓ | ✓ | ✓ | ✓ | ✓ |
| 0 Original, non-replaceable ocular | ✓ | ✓ | X | X | X |
| 2C-1.2 UV with good colour recognition, clear lens | ✓ | ✓ | X | X | X |

BCGS1 ANSI/ISEA Z87.1 rating: +Impact resistance against flying fragments, objects, large chips, particles, sand, dirt, etc.

STERILISATION:

| | BCGS1 | BVGS | BCAG | BCAP | BCAH |
|---------------|-------|------|------|------|------|
| Sterile EO | ✓ | X | X | X | X |
| Sterile Gamma | X | ✓ | X | X | X |
| Autoclavable | X | X | ✓ | ✓ | ✓ |

RECYCLABLE COMPONENTS:

| | BCGS1 | BVGS | BCAG | BCAP | BCAH |
|---|---|--|--|---|-----------------------------------|
|  Valve | Valve | Valve Pouch | Pouch | Pouch | Pouch |
|  Frame | Frame | Frame | - | - | - |
|  - | - | - | Side adjusters | Side adjusters | Vents, name plate and locking pin |
|  Polycarbonate ocular Silicone strap, PE/PET pouch | Polycarbonate ocular, Silicone strap | SEBS thermoplastic frame, Polycarbonate ocular, Silicone strap | SEBS thermoplastic frame, Polycarbonate ocular, and clip Silicone strap | Polycarbonate ocular and clip, Silicone frame and strap | |

EXPLANATION OF SYMBOLS AND PICTOGRAMS USED ON GOGGLE PACKAGING: **(01)** Please read the Instructions for Use, prior to using the goggles, or contact Nitritex for more information. **(02) CE** – Complies with the requirements of the EU PPE Regulation, (EU) 2016/425. **(03) UKCA** – Product is compliant and certified to the requirements to Regulation 2016/425 on personal protective equipment, as amended to apply in GB. **(04) SINGLE USE ONLY** – DO NOT REUSE. **(05)** Do not use if package is damaged. **WARNINGS!** - Eye-protectors against high speed particles worn over standard ophthalmic spectacles may transmit impacts, thus creating a hazard to the wearer. - If protection against high speed particles at extremes of temperature is required then the selected eye-protector should be marked with the letter T immediately after the impact letter, i.e. FT, BT or AT. If the impact letter is not followed by the letter T then the eye protector shall only be used against high speed particles at room temperature. - Materials that come into contact with the wearer's skin could cause allergic reactions to susceptible individuals. - Scratched or damaged oculars should be replaced (BCAG, BCAA and BCAP only). - To maintain the safety and effectiveness of the goggles they must be stored and transported in their original packaging. - BCAG, BCAA and BCAP, remove goggles from bag prior to autoclaving. **STORAGE** - Keep away from direct sunlight; store in a dry place and keep in the original packaging. Keep away from ozone sources. If products are properly stored, as indicated, they won't lose their performances or change characteristics significantly. If products could be affected by ageing or storage, the expiry date is mentioned on the packaging materials. **CLEANING BCAP, BCAG & BCAA GOGGLES** - Instructions for reprocessing BCAP, BCAG and BCAA goggles can be downloaded from www.ansell.com/life-sciences. Use only de-ionised or distilled water when autoclaving, autoclave at 121°C (250°F) for 30 minutes only. BCGS1 and BVGS are single use and must not be reprocessed.

- *original lens and original frames must be replaced only by the same components (replacement part codes can be found in the replacement part table)*
- *single components (frame and/or lens) must be mounted only with lens or frame of original complete product (replacement part codes can be found in the replacement part table)*
- *alone components (frame and/or lens) don't give protection*

REPLACEMENT PARTS:

| | BCGS1 | BVGS | BCAG | BCAP | BCAH |
|---------------|--------------|-------------|-------------|-------------|-------------|
| Ocular | X | X | BCAG-SL | BCAP-SL | BCAH-SL |
| Strap | X | X | BCAG-SS | BCAP-SS | BCAH-SS |

To change the ocular or strap on a goggle, simply unclip the original part and clip the replacement into its place. For ocular replacement ensure that the edge of the ocular is fully enclosed by the frame and any clips are replaced securely.

Declaración de conformidad UE disponible para su descarga en - www.ansell.com/regulatory | Organismo notificador de la UE: INTERTEK Italia S.p.A. Via Guido Miglioli 2/A, 20063 Cernusco sul Naviglio - Milano (MI), Italia. Organismo notificado N° 2575 | Organismo aprobado UKCA: ITS Testing Services (UK) Ltd, Centre Court, Meridian Business Park, Leicester, Reino Unido. Organismo acreditado N° AB0362.

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